

Custom RNA Oligonucleotides

Innovation • Value • Discovery

(V1/2015-12-14)

Bioneer Corporation

8-11 Munpyeongseo-ro, Daedeok-gu, Daejeon, 34302 Republic of Korea Tel: +82-42-930-8777 (Korea:1588-9788) Fax: +82-42-930-8688 E-mail: sales@bioneer.com Bioneer Inc.

1301 Marina Village PKWY, Suite 110 Alameda, CA 94501, USA Toll Free: +1-877-264-4300 Fax:+1-510-865-0350 E-mail: order.usa@bioneer.us.com Bioneer R&D Center

Korea Bio Park BLDG #B-702, 700 Daewangpangyo-ro Bundang-gu, Seongnam-si Gyeonggi-do, 13488 Republic of Korea Tel: +82-31-628-0500 Fax: +82-31-628-0555

Protocol for siRNA annealing

1. General Handling instructions:

The RNA-Oligonucleotides delivered by Bioneer are deprotected and purified. Although our Oligonucleotides are RNAse free, RNA is highly susceptible to degradation by exogenous RNAses introduced during handling. Therefore it is essential that all handling steps are conducted under sterile, RNAse free conditions. RNA oligonucleotides should not be handled with ungloved hands. RNAse free reagents, barrier pipette tips and tubes should be used.

Dry RNA oligonucleotides can be safely stored at -20°C for up to 6 months.

2. Annealing of siRNA:

Dissolve RNA oligonucleotides at a convenient concentration, e.g. 100 μ M, in RNAse free water. This solution should be stored at -20° C.

- Dilute each RNA oligonucleotide using annealing buffer to a final concentration of 50 μM.
- Combine 30 μl of each RNA oligonucleotide solution and 15 μl of annealing buffer. Final volume is 75 μl, final concentration of siRNA duplex is 20 μM.
- Incubate the solution for 1 minute at 90°C and cool slowly down afterwards to room temperature (over a period of ca. 45 min). Store at 4°C or on ice until ready to use.
- Annealed siRNA can be safely stored frozen at -20°C. Do not freeze-thaw more than 5 times.

3. Annealing buffer ingredients:

Annealing buffer concentration is: 30 mM HEPES-KOH pH 7.4, 100 mM KCl, 2 mM MgCl2, 50 mM NH4Ac.

BQ-042-101-04

1 www.bioneer.com

Revision: 4 (2015-11-06)